

## Protocol for preparation of hTEC/SVTERT24-B medium

page 1 of 1

Version: May 2021

Evercyte Ord. No.: CLHT-010-0024-B

---

Designation: hTEC/SVTERT24-B, human thymic epithelial cells

---

Growth medium: OptiPRO™ SFM (Gibco) supplemented with GlutaMAX™-I and G418:

Final components:

OptiPRO™ SFM (Gibco, Cat# 12309-019)

2 mM GlutaMAX™-I (Gibco, Cat# 35050-038, ready-to-use, stored at RT)

50 µg/ml G418 (InvivoGen, Cat# ant-gn5, 100 mg/ml stock solution, ready-to-use, stored at -20°C)

- take one bottle of OptiPRO™ SFM medium (500 ml) and discard 5 ml
- add 5 ml of GlutaMAX™-I to the remaining 495 ml OptiPRO™ SFM medium
- add 250 µl of G418 stock solution
  
- mix properly
- store at 4°C for a maximum of 4 weeks
- temper the medium to RT (not 37°C) before use