

Protocol for preparation of IMDM-FBS medium

page 1 of 1

Version: February 2025

Evercyte Ord. No.:	-
Designation:	TB/SVTERT350, human trophoblasts
Medium components and preparation:	<p>IMDM-FBS medium contains the herein described components and is prepared as described in the following:</p> <p>IMDM (Gibco, Cat# 12440053, stored at 4°C) 10% FBS (PAN-Biotech, Cat# P30-3031, stored at 4°C)</p> <ul style="list-style-type: none">- take one bottle (500 ml) of DMEM/F12 (1:1) and discard 50 ml- add 50 ml FBS (ready-to-use)- mix properly- store at 4°C for a maximum of 4 weeks- temper the medium to room temperature (not 37°C) before use
Related products:	<ul style="list-style-type: none">- TS medium (see <i>Protocol preparation of TS medium</i>)- TrophoUp medium (Evercyte, Cat# MHT-070), ready-to-use medium- TB/SVTERT350, human trophoblasts (Evercyte, Cat# CLHT-070-0350)